
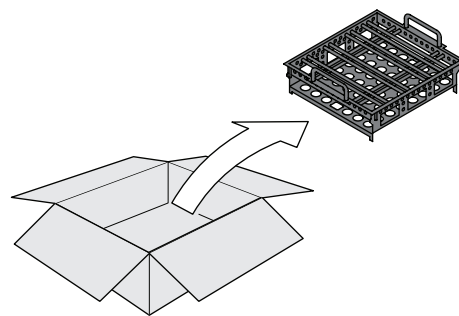
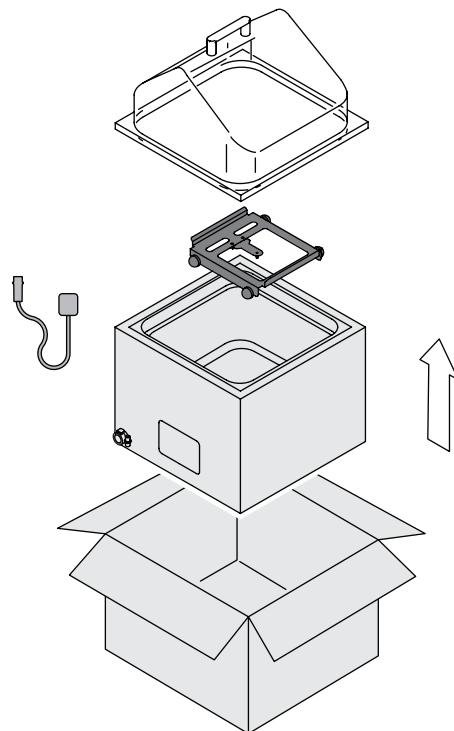
2


1

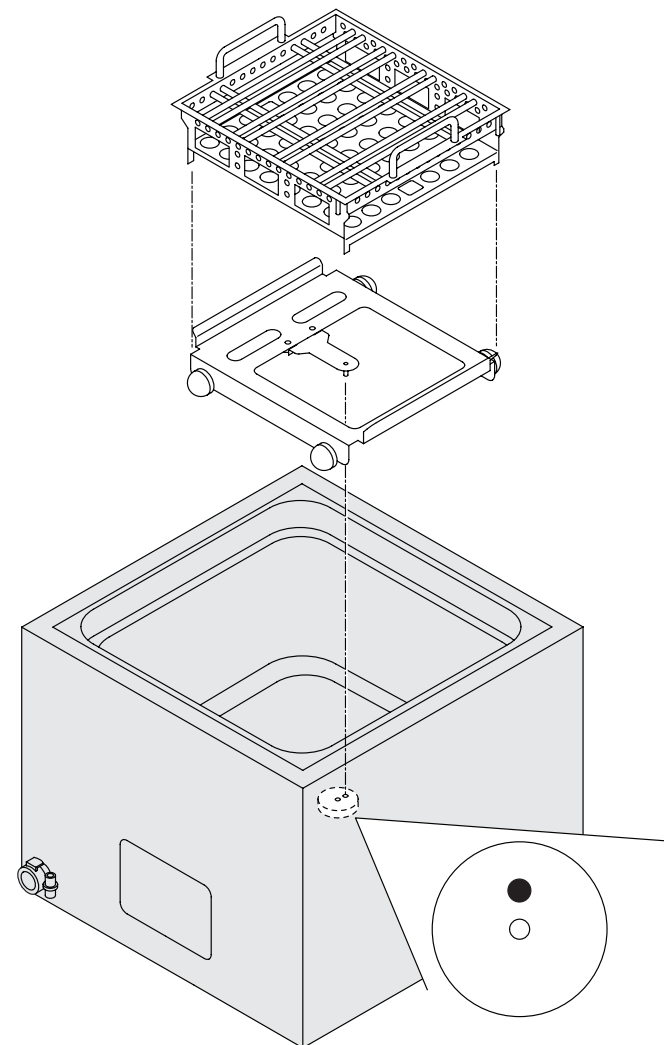
3


4



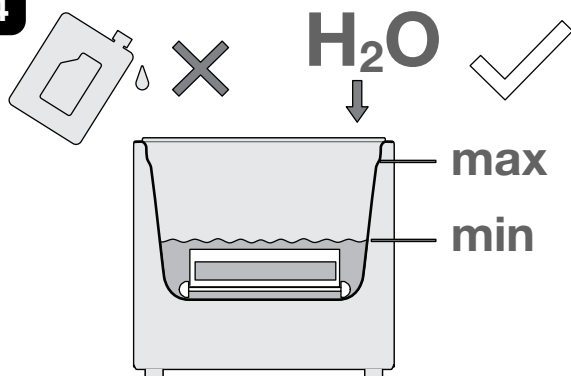
5



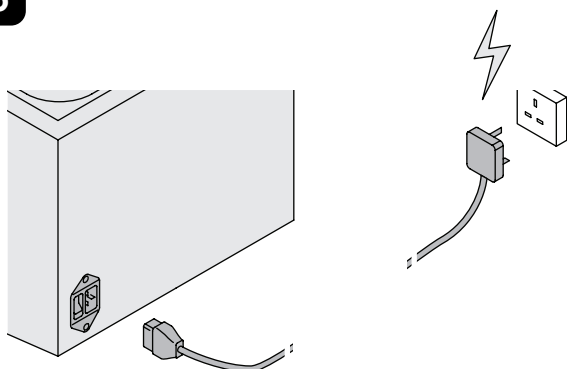
6

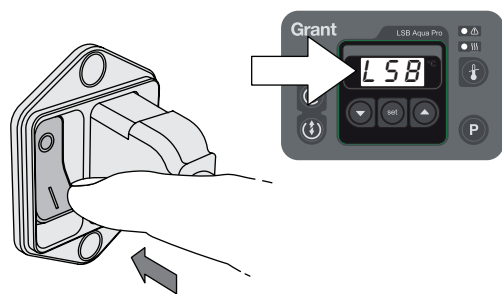
7

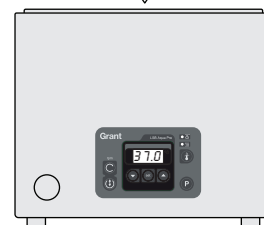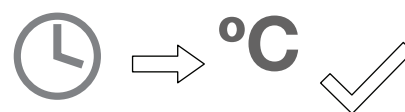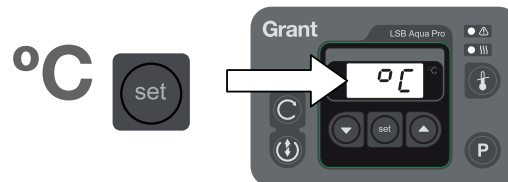
8

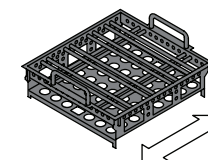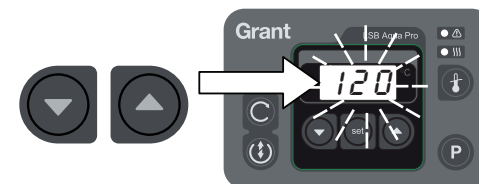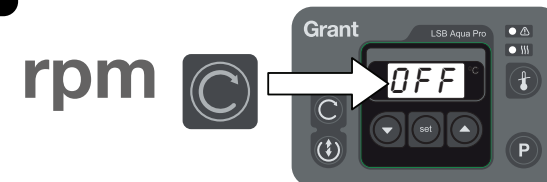stop

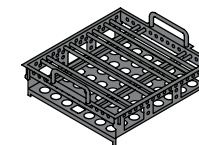
start

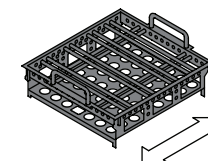
i
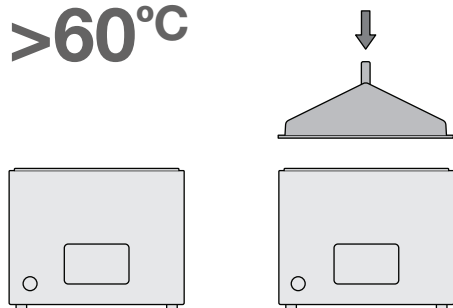
>60°C

i

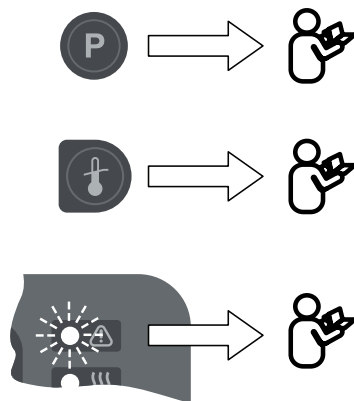Specification

Specification

Mains supply: 220 – 240V @ 50/60Hz
110 – 120V @ 50/60Hz

Earth connection required

Ambient operating temp: + 5°C to 40°C

Ambient storage temp: – 20°C to + 60°C

Current Consumption



Model	220 – 240V	110 – 120V
LSB 12	3.6A	–
LSB 12 AUS	3.6A	–
LSB 12 US	–	6.5A
LSB 18	6.2A	–
LSB 18 AUS	6.2A	–
LSB 18 US	–	9.0A

CSA mark applies to 120V models only


i
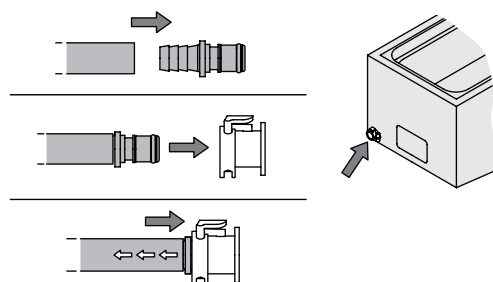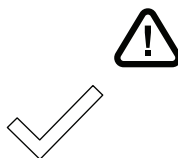
<60°C

i

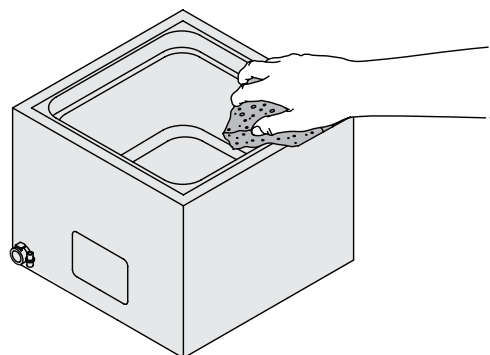H₂O ✓

Warranty & Support

3 years

When used in accordance with the user instructions this product is guaranteed for 3 years against faulty materials and workmanship.

Visit our website to register your warranty online
www.grantinstruments.com/support/warranty-registration/

Online resources
www.grantinstruments.com

Grant Instruments (Cambridge) Ltd
Shepreth
Cambridgeshire
SG8 6GB
UK

tel: +44 (0)1763 260811 fax: +44 (0)1763 262410

email: salesdesk@grantinstruments.com

Safety information

EN Original instructions

This product is a general purpose thermostatically controlled water bath designed for indoor laboratory use by or supervised by a professional user.

Read the whole of the instructions and the full Operating manual which is available online. Safety may be impaired if they are not followed.

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Surfaces and water can be hot during and after use.

This bath is only intended for use with water. Use of other fluids may invalidate warranty and present a risk of fire or explosion.

Place on a stable flat surface to reduce the risk of accidental spillage.

No user serviceable parts. Risk of electric shock after disassembly or operation with covers removed.

Not for use in environments with a risk of flammable or explosive gases.

To be operated within the limits listed below. Only use the mains cord provided. Ensure that the mains plug and the switch are easily accessible.

Use only with a mains supply which includes a ground connection.

A clearance of >10cm around the bath is required to ensure adequate air flow.

Take care when operating after a power interruption. Shaking will restart at the previously set speed.

Take care when operating with vessels in the tray. Accidental water spillage can occur when shaking is started. Reduce the water level as required.

If a potentially hazardous liquid is spilt onto the equipment, disconnect it from the power supply and have it checked by a competent person. It is the user's responsibility to carry out appropriate decontamination if hazardous material is spilt on the equipment.

There are accessible moving parts in the bath. Observe safe working practises when operating a shaking bath. Users should not interfere with moving parts when in use and prevent accidental contact from items such as fingers, hair or clothing.

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Restricting or interfering with the tray motion in the bath may reduce performance and reduce reliability.

To preserve your water bath in peak condition consult the extra guidance listed in the Operating manual. Failure to do so may affect your warranty.

Consult online resources for additional important product information.

Retain the product packaging for future storage or shipment.

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